



3 x ϕ 2.05 ∇ 6.35
M2.5 - 6H ∇ 5
 ∇ ϕ 2.55 X 90°, Near Side
EQUALLY SPACED ON 45 PCD

INVERTED (INV) MODEL IS DESIGNED TO WORK WITH THE OBJECTIVE ABOVE THE STAGE SO GRAVITY ACT IN COMPRESSION ON THE PIEZO OFTEN ON AN INVERTED MICROSCOPE.

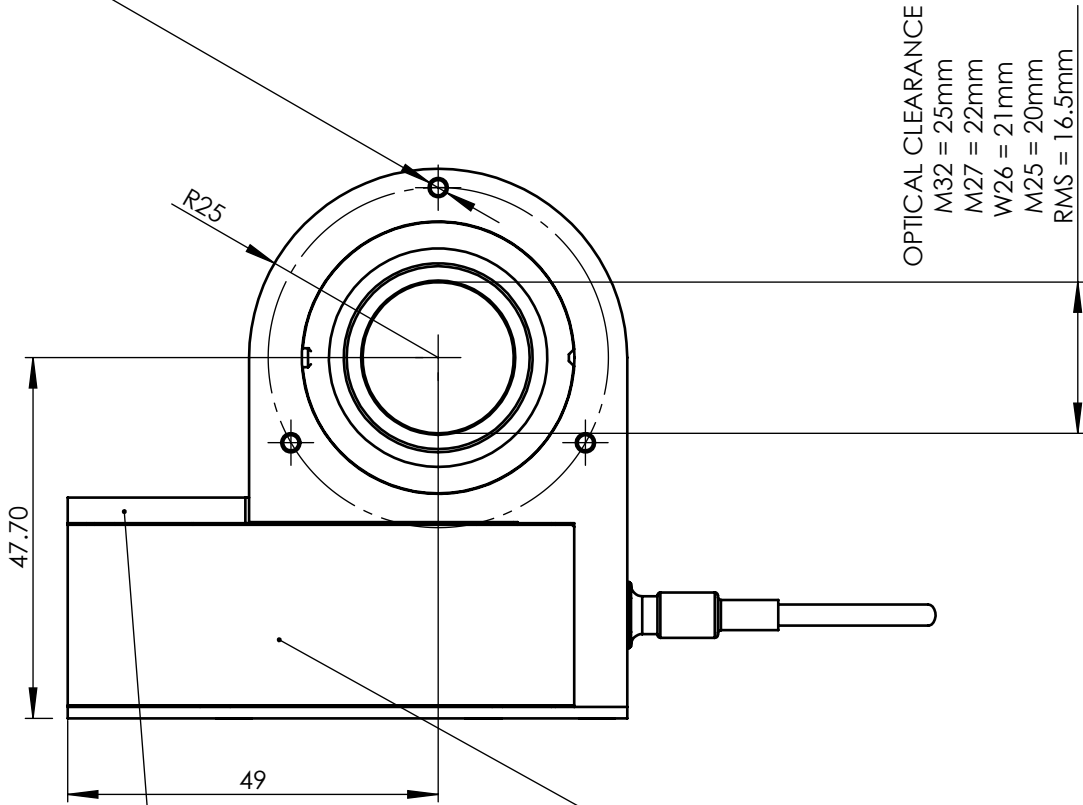


WHEN MOUNTING DIRECT ON THE 3 x M2.5 LEAVE >0.5 CLEARENCE ABOVE THIS FACE


WHEN MOUNTING DIRECT ON THE 3 x M2.5 LEAVE >0.5 CLEARENCE ABOVE THIS FACE

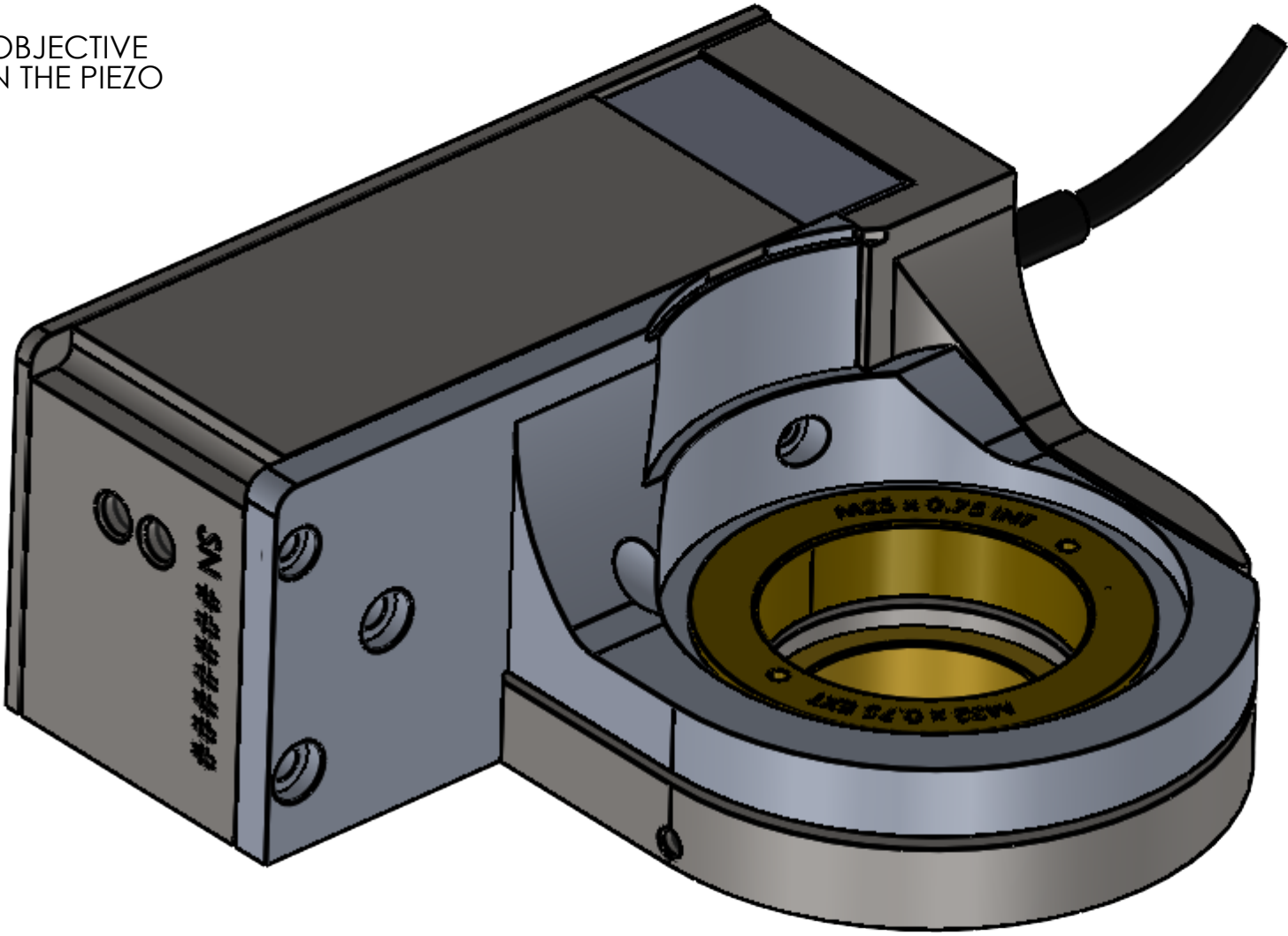

ADAPTERS AND ACCESSORIES
(SOLD SEPARATLY NOT SHOWN TO SCALE)

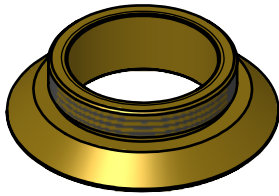


MICROSCOPE THREAD ADAPTER. SELECT APPROPRIATE THREAD TO MATCH MICROSCOPE RMS, M25, W26, M27 OR M32

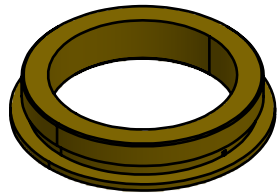

OBJECTIVE THREAD ADAPTER. SELECT APPROPRIATE THREAD TO MATCH OBJECTIVE RMS, M25, W26 OR M27. NOT REQUIRED FOR M32 SYSTEMS.




PARAFOCAL SPACERS FOR MOVING OTHER OBJECTIVES IN MICROSCOPE NOSE TO THE SAME FOCAL HEIGHT RMS, M25, W26, M27 or M32. 13mm LONG EXCEPT M32 WHICH IS 12mm.

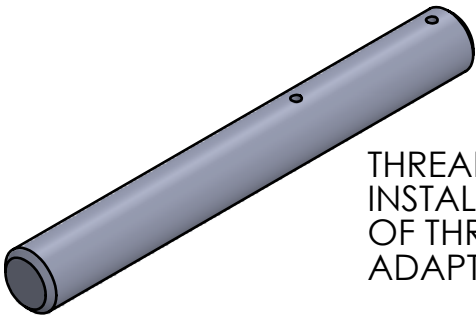


THREAD ADAPTER KEY CAN AID INSTALLATION AND REMOVAL OF THREAD ADAPTERS LIGHTLY NIP UP ADAPTERS ON INSTALLATION.

SELECT MICROSCOPE THREAD ADAPTER WITH THREAD TO SUIT MICROSCOPE

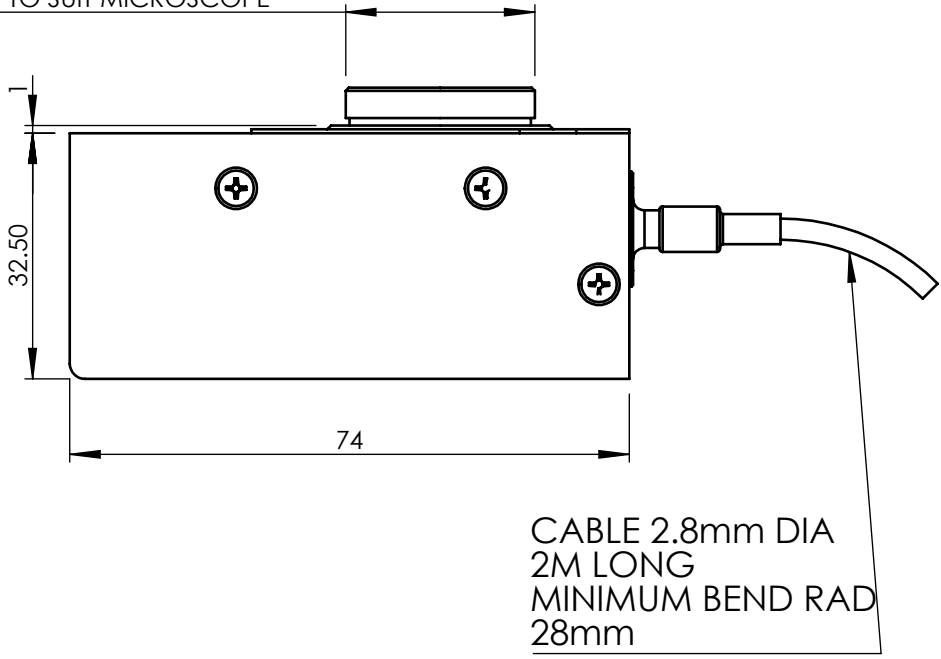


CABLE 2.8mm DIA
2M LONG
MINIMUM BEND RAD
28mm

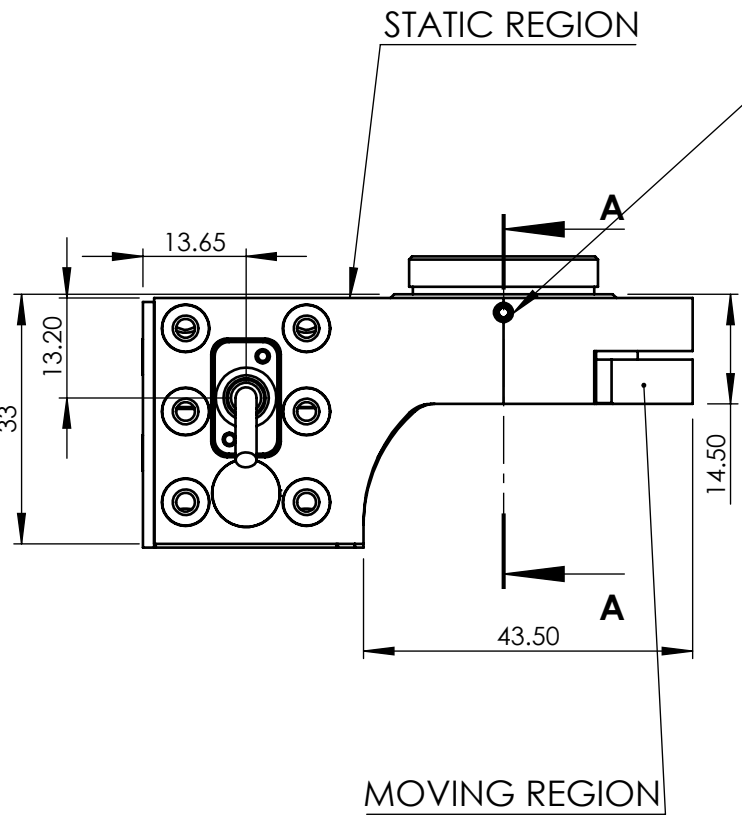

STATIC REGION

MOVING REGION

TWO MICROSCOPE ADAPTER CLAMPING GRUB SCREWS USE 1.3mm (0.050") HEX KEY. USE ONLY THE SCREWS PROVIDED.

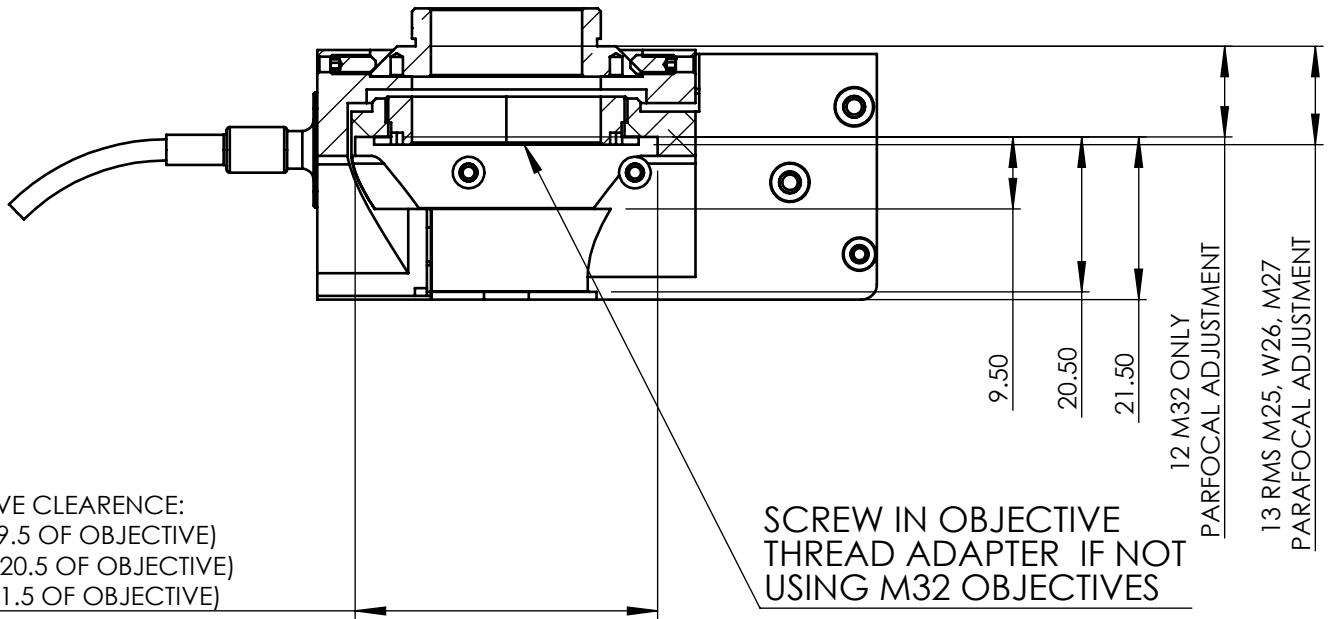

OBJECTIVE CLEARENCE:
 ϕ 40 (TOP 9.5 OF OBJECTIVE)
 ϕ 42.8 (TOP 20.5 OF OBJECTIVE)
 ϕ 44 (TOP 21.5 OF OBJECTIVE)

SCREW IN OBJECTIVE THREAD ADAPTER IF NOT USING M32 OBJECTIVES

SECTION A-A

TITLE	OP800-INV INSTALLATION DRAWING	APRVD.	S.R.	3RD ANGLE  DO NOT SCALE IF IN DOUBT ASK
DRG. No.	QG-0494-IN-100004	DRAWN	JC	
MATL.	STAINLESS, BRASS, ALUMINIUM	SCALE	1:1	
FINISH	CLEAN	DATE	02.01.25	
PRIOR Scientific Instruments Ltd. Fulbourn, Cambridge.				TOLERANCE UNLESS OTHERWISE STATED $\pm$ 0.1mm BREAK ALL SHARP EDGES

ECN	REV	CREATED	APP	DATE	DETAILS
1st Rel	B	JC	S.R.	06.01.25	1st Release